

Agricultural chemical Analysis by HPLC

1. Analytical Condition

YL9100 HPLC

Mobile Phase : Water, ACN Gradient

Flow Rate : 1.5ml/min

Column Oven : 40°C

Detector : UV/Vis 254nm

Column : Kinetex C18 (4.6*100mm, 2.6um)

Injection Volume : 10ul

<Gradient Program>

Gradient Program	Water	ACN
Initial	90	10
15.00	10	90
16.00	10	90
16.10	90	10
25.00	90	10



Agricultural chemical Analysis by HPLC

2. Chromatogram

